

# ZD-8800

## Horizontal Receptacle Shaker

### Operational Manual

Omxie Corp.  
<http://www.smtmax.com>



## **1. Technical Data**

Shaking mode: alternating motion  
Shaking speed : 10-280 (rpm)  
Display: digital  
Unit size: 440×390×230(mm)  
Plat-form size: 400×360(mm)  
Clamping fixture: spring net  
Power: AC110V 50-60HZ  
Net weight: 25kg

## **2. Installation**

- The shaker should be placed on a flat platform.
- There should be no other objects around it within 10-20cm.
- The surrounding temperature should be between -10°C—40°C.
- The shaker should not be used in wet or high-temperature environment.
- Make sure good power and ground connection.

## **3. Operations**

1. Press the "Power" button. The power light and LED's should be on to indicate that the shaker is ready.

2. Adjust the shaking speed: turn the "speed" knob clockwise to make it faster; turn it counterclock to make it slower. The display window changes accordingly to indicate the shaker's current speed.

3. Adjust height of spring net clamping. Turn the black handwheel on the side of the net to set it to one of the two available levels.

Contact:

Omxie Corp.  
125 Business Center Dr Unit G  
Corona, CA 92880  
Toll Free: 1-951-543-4543  
<http://www.omxie.com>