

# ZD-9550 ROCKING SHAKER Manual


Omnie Corp.  
<http://www.smtmax.com>



## 1. Technical Specifications

Shaking speed: 20-65 rpm  
Display: LED  
Swinging range:  $\leq 20^\circ$   
Platform size: 10.6"×10.6"  
Max load capacity: 10kg  
Dimensions: 10.4"×10.4"×6"  
Supply power: 110V AC, 40W, 50-60HZ

## 2. Front Panel

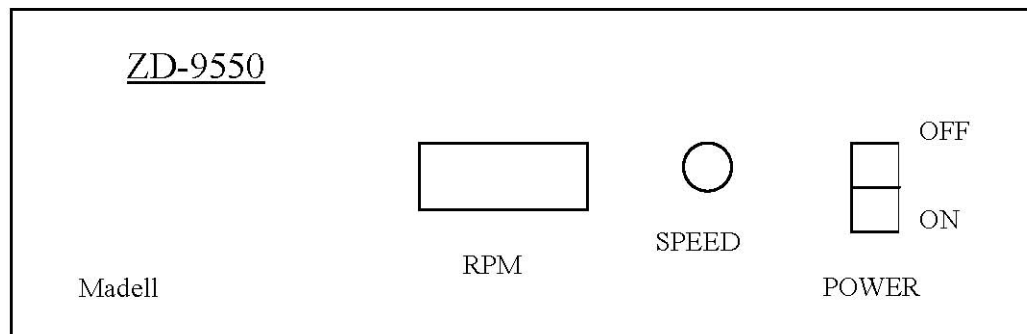

## 3. Operations

1. Connect the power supply; turn on the "Power" switch. The power LED is on to show that the shaker is ready.
2. Adjust the shaking speed: turn the "speed" knob clock-wise makes the shaker faster, counterclock wise makes it slower. The "RPM" display changes accordingly to reflect the new speed.
3. Adjust the swinging angle: this function is located in the middle of the shaker's back panel. **Before adjusting, turn off the power supply and make sure that the shaker has stopped completely. Otherwise, it may cause serious finger harms.** Loosen the roller screw counterclock wise, turn the angle adjustment handle to the proper position, and tighten the screw.

Contact:

Omxie Corp  
125 Business Center Dr Unit G  
Corona, CA 92880  
USA  
1-951-543-4543  
<http://www.smtmax.com>